

---

## ABOUT US

---

Molecular Toxicology, Inc. is the leading manufacturer of products used in the Salmonella and E. coli WP2 mutagenicity tests. Moltox minimal glucose agar plates, top agars, Salmonella and E. coli tester strains, frozen and lyophilized S9, MUTAZYME™, NADPH-regenerating systems and positive control chemicals are distributed worldwide. Moltox has developed microtiter plate format fluctuation tests consistent with OECD guidelines including Moltox® FT™ tests and distributes the BioReliance Ames II™ test kit.

# MOLTOX®

Molecular Toxicology, Inc.

---

### MOLTOX Distributors

---

#### EUROPE

Trinova Biochem GmbH  
[www.trinova.de](http://www.trinova.de)

#### INDIA

Krishgen Biosystems  
[www.krishgen.com](http://www.krishgen.com)

#### PUERTO RICO

Instrumed  
[www.instrumed.net](http://www.instrumed.net)

#### SOUTH KOREA

Woo Jung BSC, Inc.  
[www.woojungbsc.co.kr](http://www.woojungbsc.co.kr)

#### CHINA

Shanghai Bioplus Biotech Co., Ltd.  
[www.bioplus-biotech.com](http://www.bioplus-biotech.com)

#### BRAZIL

Interlab  
[www.interlabdis.com.br](http://www.interlabdis.com.br)

#### NORTH AMERICA

Brand-Nu Laboratories  
[www.brandnu.com](http://www.brandnu.com)

Fisher Scientific  
[www.fishersci.com](http://www.fishersci.com)

VWR Scientific  
[www.vwrsp.com](http://www.vwrsp.com)

---

### Contact Us

---

Molecular Toxicology, Inc.  
157 Industrial Park Drive  
PO Box 1189  
Boone, NC 28607 USA

Tel: 828.264.9099  
Toll Free: 800.536.7232  
Fax: 828.264.0103  
Email: [Sales@MOLTOX.com](mailto:Sales@MOLTOX.com)

[www.MOLTOX.com](http://www.MOLTOX.com)


# MOLTOX®

Molecular Toxicology, Inc.

---

## MOLTOX® FT™ TA98/TA100 Mutagenicity Assay Kit #31-300

---

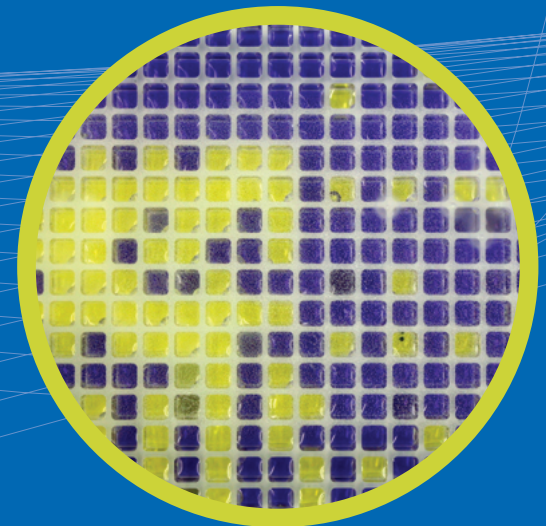

# MOLTOX<sup>®</sup> FT<sup>™</sup> TA98/TA100 Mutagenicity Assay Kit


## Kit Components

Kit #31-300 includes:

PART # DESCRIPTION

71-098.2L	TA98 inocula (2 discs/vial)
71-100.2L	TA100 inocula (2 discs/vial)
26-712.05	Moltox <sup>®</sup> FT <sup>™</sup> Growth Medium (50mL)
26-710.05	Moltox <sup>®</sup> FT <sup>™</sup> Exposure Medium (50mL)
26-711.3	Moltox <sup>®</sup> FT <sup>™</sup> Reversion Indicator Medium (300mL)
11-401.3	MUTAZYME <sup>™</sup> (30% s9 Mix, 3.25mL)
60-157	2-Aminoanthracene (100µg/vial)
60-159	4-Nitroquinoline-N-oxide (50µg/vial)
60-161	2-Nitrofluorene (50µg/vial)

User Instructions Included

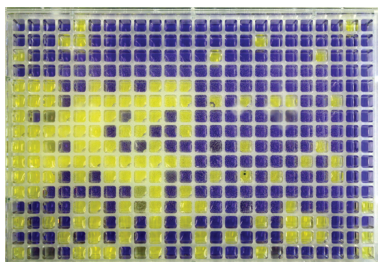

384-Well  
Microtiter Plate

## Basis of the Test

Moltox<sup>®</sup> FT<sup>™</sup> microplate fluctuation tests estimate the mutagenic potential of a test material by measuring its ability to induce reversion of *his*<sup>-</sup> *Salmonella typhimurium* or *trp*<sup>-</sup> *Escherichia coli* auxotrophs to their respective prototrophic conditions. The bacterial strains used in the assays are identical to those used in conventional plate incorporation assays (Mortelmans, K and E Zeiger, *Mutat res*, 455:29-60, 2000 and Mortelmans, K and E Riccio, *Mutat res*, 455:61-69, 2000). The experimental design used in the Moltox<sup>®</sup> FT<sup>™</sup> tests is based upon D. Gatehouse's adaptation of the design reported by S. Luria and M. Delbruck in 1943 (see, *Mutat res*, 53:289-296, 1978; Genetics, 28:491-511, 1943). *Trp*<sup>-</sup> or *his*<sup>-</sup> target cell populations are treated in 24-well plates using media containing limiting L-tryptophan or L-histidine. After the treatments, the cells are transferred to 48-well sectors of 384-well microtiter dishes in a L-tryptophan-L-histidine-free medium containing a pH indicator. After 48 hours incubation, cells able to grow in the absence of L-histidine or L-tryptophan (mutant cells) will have proliferated resulting in media acidification and the appearance of yellow wells. These wells are counted and the number for each treatment condition compared to their negative control (Gilbert, RI, *Mutat res*, 74:283-289, 1980).



## Advantages of the Test

- Fluctuation tests in microtiter plate format exhibit increased sensitivity and may be automated.
- By combining certain strains, costs can be reduced and test material conserved.
- Assays are conducted using activation and non-activation conditions similar to those employed in the traditional Ames test.
- Colorimetric endpoint is unambiguous and results are easily submitted to statistical analysis.
- Complies with OECD 471 Guideline for Testing of Chemicals (Bacterial Reverse Mutation Test).

