


Culturing of MOLTOX[®] Lyophilized Discs


Day 0 (end of day)

Day 1


VIA FLASK

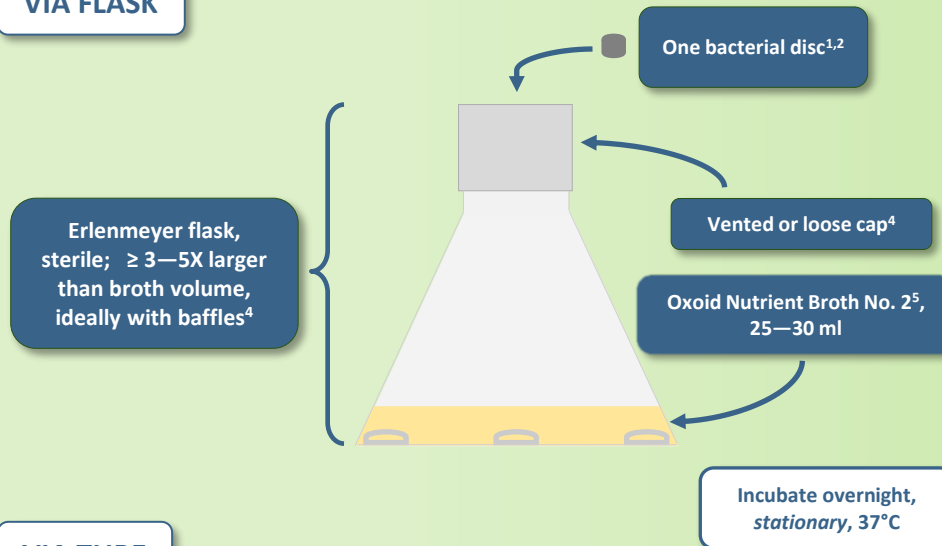

VIA TUBE

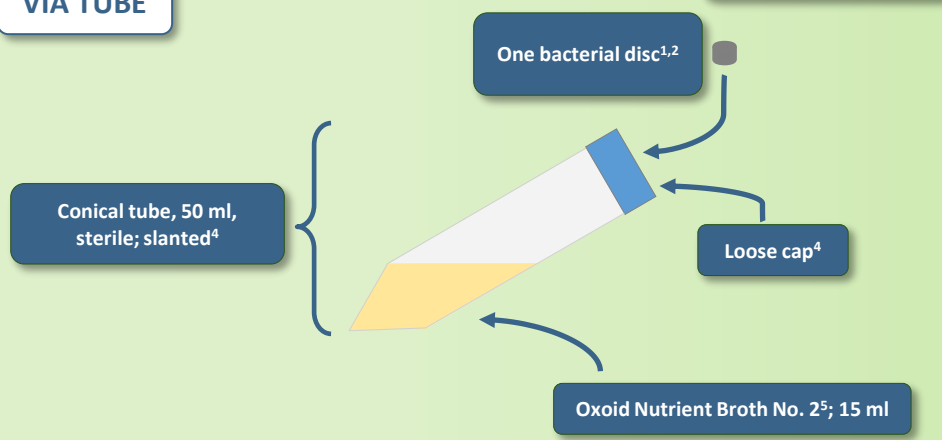

AM (upon arrival)

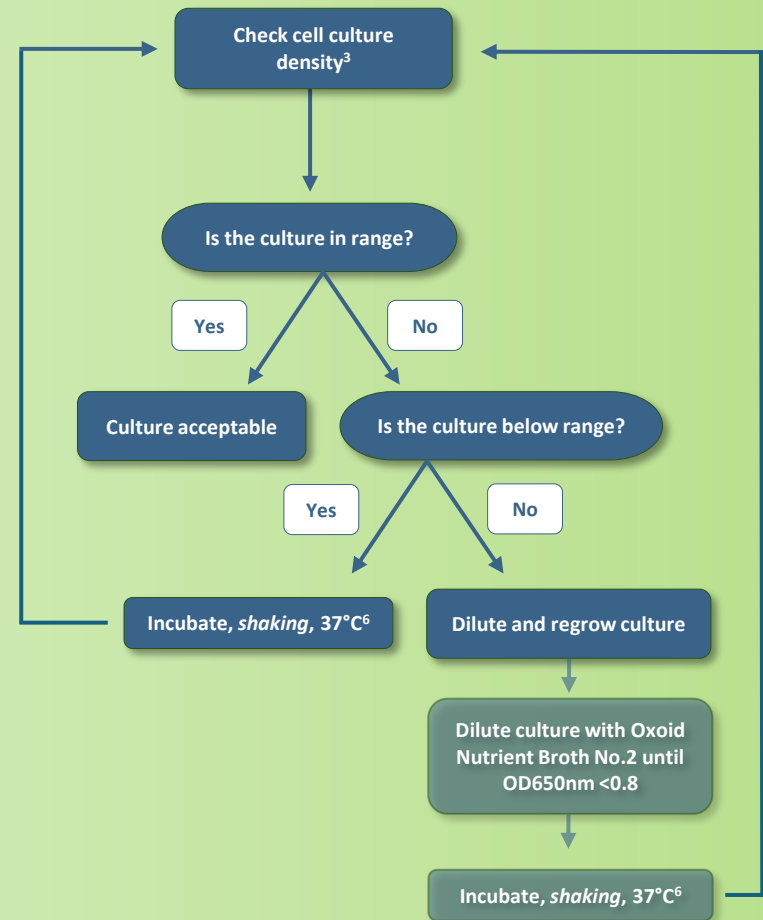

¹ Also reference MOLTOX[®] Technical Bulletin *Lyophilized Culture Discs*.

² Allow vial to come to room temperature prior to opening. Seal tightly after use. Store refrigerated.

³ OD650nm; 1.0–1.4. Verification via an internal growth curve study is recommended.

⁴ Increased surface area, loose/vented caps, and baffles increase aeration leading to better growth.

⁵ Oxoid Nutrient Broth No. 2 is the **only** acceptable growth medium.

⁶ Incubation/shaking time dependent on initial OD reading. Strains can double every 20 minutes.